



## Chiral Analysis of Flurbiprofen using a Circular Dichroism Detector

### Introduction

Flurbiprofen is a non-steroidal anti-inflammatory drug and used as a topical medical patch or an internal medicine. The S enantiomer of flurbiprofen, accounts for the majority of the pharmacological activity.

The Circular Dichroism detector is well known for its high accuracy and selectivity for chiral analysis of compounds that have chromophores close to the chiral center. Both CD and UV signal are measured simultaneously and spectral scanning can be used to optimize the wavelength of interest, which is important in the measurement of chiral compounds. Not only can the CD detector help differentiate between two enantiomers, but it can also distinguish between achiral impurities and enantiomers of interest.

**Keyword:** Flurbiprofen, enantiomer, CD detector, CHIRALPAK AD-RH


### Experimental

#### [Equipment]

Pump: PU-4185  
 Pump option: Degasser unit  
 Autosampler: AS-4050  
 Column oven: CO-4060  
 Detector: CD-4095

#### [Conditions]



Column: CHIRALPAK AD-RH (4.6 mm I.D. x 150 mm L, 5 µm)  
 Eluent A: 0.1% phosphoric acid aqueous solution  
 Eluent B: Acetonitrile  
 Composition: Eluent A/B (50/50)  
 Flow rate: 0.5 mL/min  
 Column temp.: 25°C  
 Wave length: 250 nm  
 Response: 1.5 sec  
 Scan speed: 10 nm/sec  
 Injection volume: 10 µL  
 Standard sample: 100 µg/mL racemic flurbiprofen in eluent A/B (50/50)

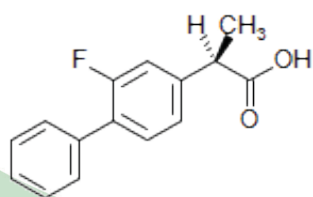

Flurbiprofen

## Results

Fig. 1 shows the chromatograms of the racemic flurbiprofen standard. Fig. 2 CD and UV spectra (measured using stopped-flow).

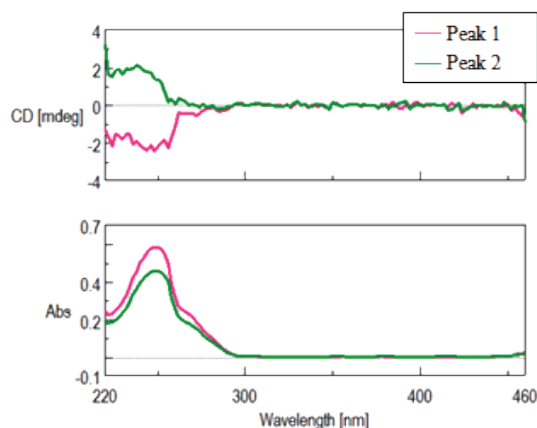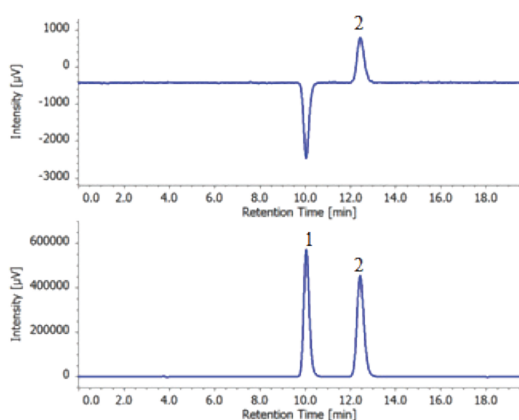

**Figure 1.** Chromatogram of Flurbiprofen standard  
Top: CD detection, Bottom: UV detection

**Figure 2.** Spectral measurement Top: CD, Bottom: UV

P/N	Description
7003-J014A	PU-4185 RHPLC Semi-micro Pump (Base Unit)
7006-H003A	DG-4000-04 4-channel Degasser Unit, for Analytical
7017-J002A	AS-4050 HPLC Autosampler
7021-J002A	CO-4060 Column Oven
7029-J002A	CD-4095 Circular Dichroism Chiral Detector
7058-J011A	BS-4000-1 Bottle stand
6688-H564A	LC-Net CG cable
7059-J012A	ChromNAV Ver.2 Chromatography Data System
7001-H401A	HPLC Start Up Kit for LC-4000
7001-H405A	Maintenance tool kit
	19724 CHIRALPAK AD-RH manufactured by Daicel